

VFC & VFAAR




DEPARTMENT OF
PUBLIC HEALTH

Identifying VFC-400 Alarms


If your digital data logger's alarm goes off, you might have an out-of-range temperature that puts vaccines at risk.

Here's how to identify an alarm



No alarm

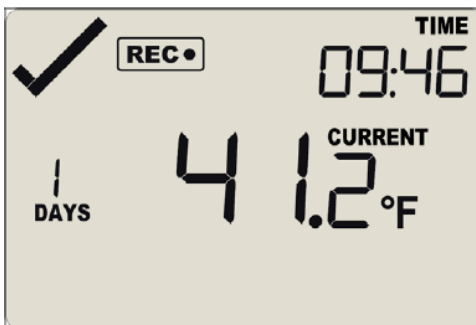

Alarm

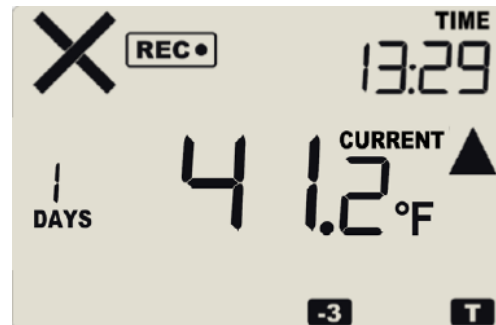

Can you spot the difference?

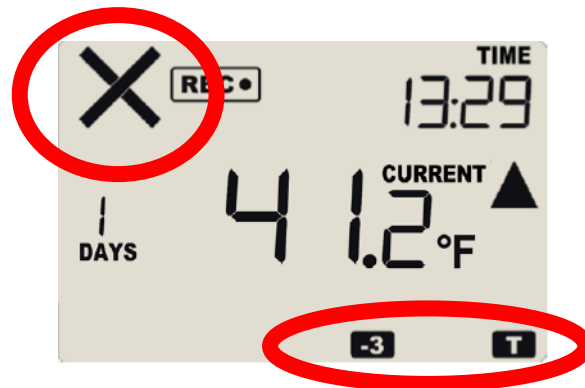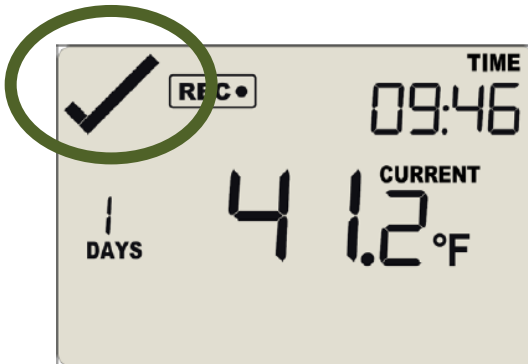

No alarm

A checkmark in the upper-left corner

No black boxes below temperature

Alarm

X in the upper-left corner

Black boxes below the temperature

New Feature: The VFC-400 includes audible alarm. The unit will beep when there is an alarm.