

DDL Stop/Start Guide

Storage & Handling

Questions about your Digital Data Logger (DDL)? Contact our Storage and Handling team at TempCheck@phila.gov or 215-685-6777

Stopping the DDL

Step One: **Stop your DDL**

- Press and hold the *Start/Clear/Stop button* until the word **STOPPED** appears on the screen (fig. 1).
- When **REC•** disappears, release the *button* immediately.
- Your DDL is now in **stopped mode** (fig.2).

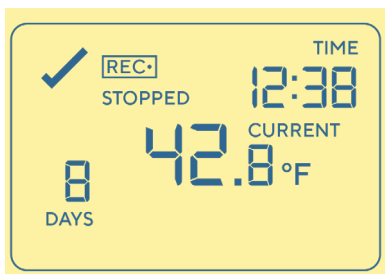


fig. 1



fig. 2

Step Two: **Disconnect the Probe**

- Unplug your DDL from its probe (fig.3).

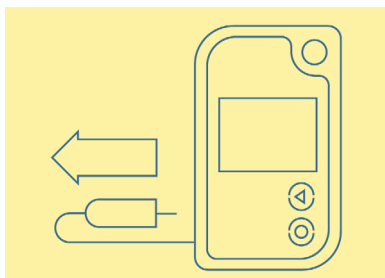


fig. 3

Step Three: **Download Data**

- Take your DDL to your computer, open the software and insert the logger in the docking station to download the data.
- Repeat the process every 30 days.

DDL Stop/Start Guide

Storage & Handling

Questions about your Digital Data Logger (DDL)? Contact our Storage and Handling team at TempCheck@phila.gov or 215-685-6777

Starting the DDL

Step One: **Connect Your DDL to the Probe**

- Reconnect the probe to your DDL (fig. 4)

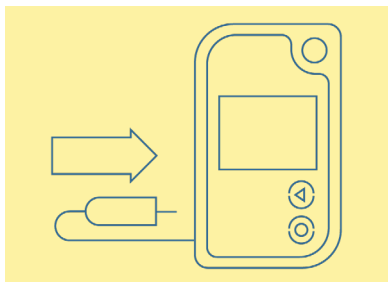


fig. 4

Step Two: **Start Your DDL**

Your DDL must be in **ready mode** to start as shown in fig. 5. If your DDL is not in ready mode, please contact TempCheck@phila.gov for troubleshooting.

- Press and hold the *Start/Clear/Stop button*. **STARTING** will appear (fig. 6).
- When **READY** disappears, **release the button immediately**.
- **STARTING** will be replaced with **REC•** and the current temperature reading will appear (fig. 7). **Your DDL is now recording temperature!**

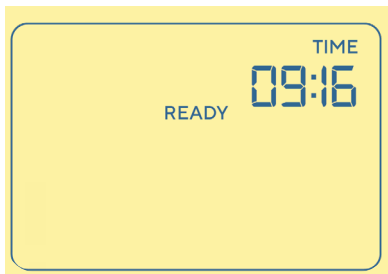


fig. 5

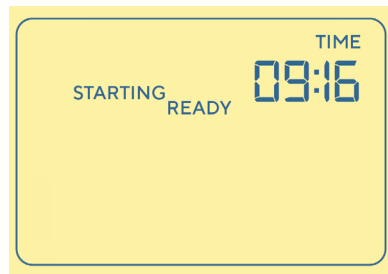


fig. 6

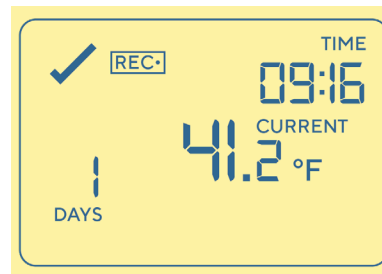


fig. 7